Preparative Module 6100

V 03/2006

The **Preparative Modul 6100 V03** is based on the V315 variant of the Assistant 6000. It is equipped with a Smartline Pump 100 and an S-6 valve drive on the front side and with an S-12 valve drive in the rear. The Module 6100 includes a 50 ml pump head, an injection valve, a fractionation and recycling valve and the preparative accessory kit. All capillary connections are designed for 1/8" capillaries.

By combining these single modules into the **Smartline Assistant 6000** unit, these modules are seamlessly integrated into a Prepline system or a Smartline tower. Operation and control of the single modules is not affected by integration into the Assistant 6000 with one common power supply.

The only exception to this is that the valve drive in the rear cannot be operated manually but only via software or via the remote control strip and event signals given by other controlling instruments. The control connectors (RS-232 and remote control strips) are located on the rear side of the Assistant 6000.

The **Preparative Module 6100** is offered in several configurations, for instance with either 6, 12, or 16 switching positions on the fractionation valve, with either 1/8" or 1/16" connectors.

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The **fractionation/recycling valve** enables the collection of up to 5, 11, or 15 fractions. Port 6, 12, and 16 are the waste positions, respectively. Optionally, two ports can be used for solvent and peak recycling.

The number of fractions can be easily expanded by cascading one or more additional valves. The position numbers are added sequentially in the order that the additional valves are connected. The last port of a valve is used to connect to the central input connector of the subsequent valve. The waste and recycling positions are always situated on the last valve.

To achieve a clear tubing arrangement, the **guidance metal sheet** should be used. It can be placed between two Smartline or Prepline instruments either with the narrow or with the broad side facing out. The feet on the left-hand side of the upper instrument are to fit into the sheet recesses.

Rubber feet on both sides of the sheet guarantee stable positioning.

The large amount of drill holes makes it possible to easily organize fractionation with collection vessels of almost any size or shape.

Exchange of fraction collection racks is made easier by not guiding the tubings into the vials. The accurate dropping of fractions into collection vessels is done by using a stainless steel adapter affixed to the vial rack.

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Technical data are subject to change without notice.

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Visit www.knauer.net for details on complete HPLC systems, HPLC columns, and osmometers.

KNAUER · ASI · Advanced Scientific Instruments
 Wissenschaftliche Gerätebau
 Dr. Ing. Herbert Knauer GmbH
 Hegauer Weg 38, 14163 Berlin, Germany
 Telephone: +49 (0)30 / 80 97 27-0
 Telefax: +49 (0)30 / 8 01 50 10
 E-Mail: info@knauer.net
 Internet: www.knauer.net

Your local distributor: