Euroline

Eurosil Bioselect

Eurosil Bioselect – special material for separation of biopolymers

Eurosil Bioselect 300 phase material was developed for the separation of proteins and peptides. By derivatisation of special pretreated large pore size silica gel Eurosil Bioselect material is available with long chain C18 modification and short chain C4 modification.

The quite close and strongly controlled pore distribution is a main quality parameter as the absence of interfere anions and heavy metal ions on the silica surface is constantly verified.

With a pore size of 300 Å the specific surface (BET) of the Eurosil Bioselect phase is calculated with 100 m²/g and the pore volume is 0.8 ml/g. The carbon content of the Eurosil Bioselect phase depends on the modification. The C18 modification has a carbon content of 7.5 % and is most suitable for separation of hydrophilic biomolecules and peptides.

The second available Eurosil Bioselect phase is characterised by a short chain C4 modification. This is preferable used for the separation of hydrophobic biomolecules and large proteins. The C4 phase has a carbon content of 2 %.

The Eurosil Bioselect 300 phase is available with particle size of 5 µm, suitable for analytical and preparative purposes.

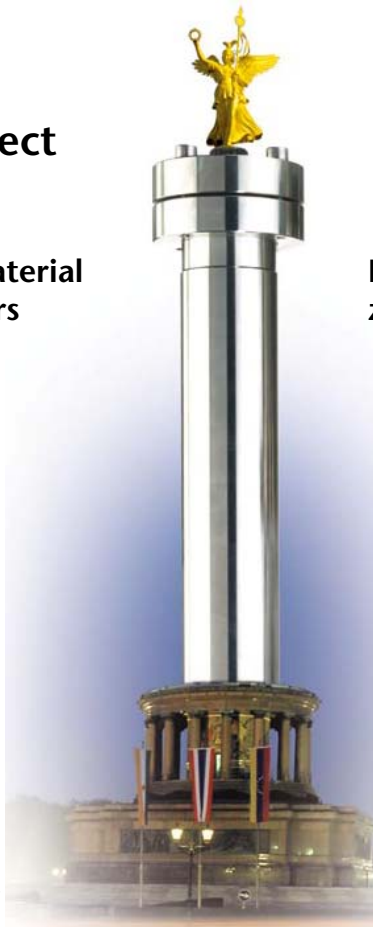
Fields of Application:

Reversed-Phase-Chromatography

- Proteins up to MW approx. 250 000 Dalton especially hydrophobic substances such e.g. membrane proteins
- Peptides especially „peptide-mapping“ after enzymatic cleavage
- Oligonucleotides

Characteristics:

- total surface coverage with C4 or C18 groups
- complete endcapping
- no non specific interaction with the silica gel matrix
- no irreversible adsorption
- maximum substance recovery
- high separation selectivity and chromatographic resolution
- very good loadability
- long life-span

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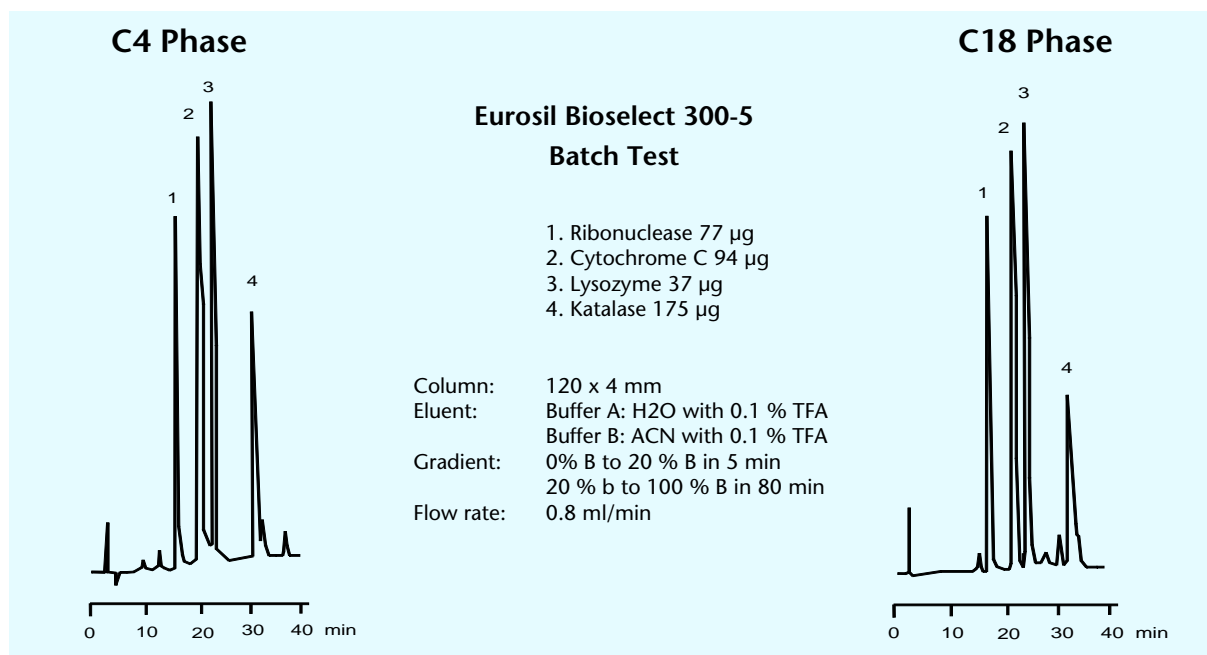
Quality Control

Every column that leaves our production plant has been individually tested and is furnished with a certificate of quality. We guarantee the quality of every HPLC column we deliver. Our quality standards with regard to peak symmetry and plate counts are extremely high to ensure consistent material and column quality.

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Other column sizes are available upon request.

Technical data are subject to change without notice.

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Visit www.knauer.net for details on complete HPLC systems, HPLC columns, and osmometers.