Smartline

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An important aspect for quality as well as for time saving concerning the analytic is the sample preparation. The Sample Preparation Unit 6300, a combination of the Smartline Assistant 6010 with the Smartline Pump 1000 (refer separate brochure) automates and minimizes the solid phase extraction process (SPE) in routine analysis and method development.

All steps of the conventional solid phase extraction can be automated with the Smartline Assistant 6010:

- Conditioning
- Equilibration
- Sample task onto the cartridge
- Elimination of matrix
- Drying of the packing material of the cartridge by a compressor
- Elution of the analyte by In-Line-Transfer to the LC System

Realise an online solid phase extraction with direct elution onto the HPLC column. As a result, the reliability of analytical data will be increased. While the HPLC part performs the chromatographic separation, the sample preparation module prepares the next sample.

Take advantage of the online solid phase extraction directly coupled with the LC System:

- Extraction with high sample throughput
- Minimized solvent consumption
- A fully automated system, thus no manual operation required by the analyst
- High reproducibility and constant flow rates
- Minimized carryover of samples due to controlled washing steps
- Fast method development of solid phase extraction process
- Simple integration into an existing LC-System by „Stand-alone“-Mode

The Smartline Assistant 6010 combined with the Smartline pump 1000 perfectly fits your needs in the fields of

- Pharmaceutical chemistry
- Diagnostics
- Forensic Analysis
- Food chemistry
- Environmental chemistry

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Setup and operation of the Smartline Assistant 6010

The Smartline Assistant 6010 includes two 6-port/3-channel-valves and one gradient mixing block, normally part of the Smartline Manager 5000. It guarantees highest gradient- and flow precision. Both valves as injection- and bypass valves are controlled by the pump.

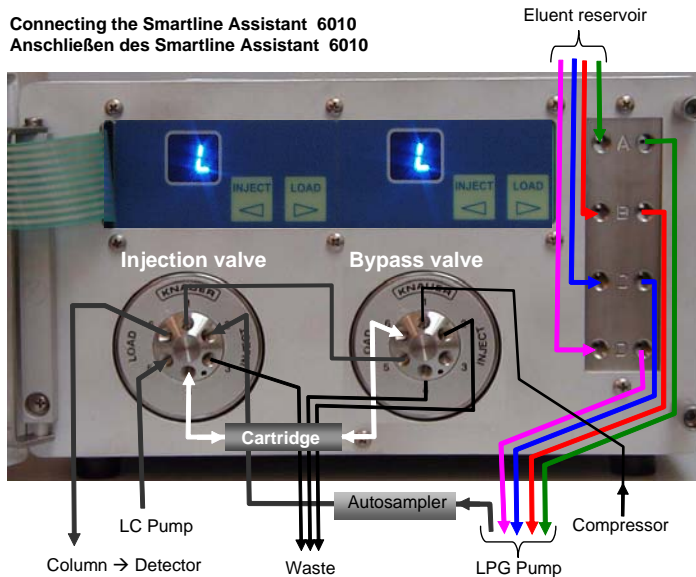
Coupling of the Smartline Assistant 6010 to a LC System

The Smartline Assistant 6010 can be combined as well with a KNAUER Smartline HPLC system as with any other HPLC system.

LC pump, HPLC column, outlet of the gradient mixing block and one side of the cartridge are connected to the injection valve. The other side of the cartridge is connected to the bypass valve. This offers the possibility to fill all capillaries with new eluent to avoid interferences mixing different eluents. Additionally the cartridge can be dried by a compressor.

The sample can be presented onto the cartridge by means with an autosampler. The autosampler is connected to the gradient mixing block and the injection valve. If higher amounts of sample should be enriched onto the cartridge as it is possible with an autosampler, the sample can be presented directly by one channel of the gradient mixing block.

Connecting the Smartline Assistant 6010
Anschließen des Smartline Assistant 6010

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Technical Data

Power supply	90-260 V, 47-63 Hz, 100 W
Dimensions	226 x 135 x 390 mm (WHD)
Weight	3.5 kg
Gradient	Number of eluents up to 4
	Range 0 – 100 %
	Graduation 1 % Steps and linear
Valves	Materials Stator: Stainless steel Rotor: Vespel
	Pressure range up to 30 MPa
	Wholes 1/16" – Connectors
	Switching time < 150 ms
	Control Short-circuit contact

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Ordering information

Order No.	Article
A5003	Sample Preparation Unit 6300 consisting of
A5003V10	Smartline Assistant 6010,
A50301	Smartline Pump 1000 with 10 ml Pump head
A5326	Eluent tray

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Technical data are subject to change without notice.

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Visit www.knauer.net for details on complete HPLC systems, HPLC columns, and osmometers.