Smartline

Smartline – HPLC by Design

Modern technology and practical design are prominent features of the Smartline series of HPLC instruments. Its thoughtful design not only conserves valuable bench space, but also allows the length of the capillaries between instruments to be kept to a minimum, resulting in a marked decrease in the dead volume of the complete HPLC system.

The display panel is integrated into the front door which hides components such as the pump head and flow cell from view. These important elements are also thereby protected against external influences. The complete HPLC system appears much more organized and attractive. The uniform Smartline design offers extended adaptability for flexible combination of individual modules.

Whether you would like to put together a system for isocratic or gradient HPLC, or an HPLC system with more than one detector, KNAUER Smartline can accommodate all of your requirements.

Our offer to you:

- ▶ You let us know which substances you would like to separate or analyze.
- ▶ We work together with you to find the right application and the corresponding column.
- ▶ We will then put together a complete HPLC system for you from our Smartline components which is best suited to your specific needs so that you only buy what you really need.

The following is a brief overview of our assortment of Smartline HPLC instruments, including autosamplers, pumps and detectors, as well as software.

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Smartline – HPLC by Design

Smartline Pump 1000

Our versatile dual-piston pump offers outstanding flow rate precision ($RSD \leq 0.1\%$) and accuracy. Its optimized mechanics provide for a long lifetime.

Up to four different eluents can be mixed in a quaternary gradient in either a low pressure (LPG) or a high pressure gradient (HPG) system. For an LPG system, the Smartline Manager 5000 is required while for an HPG system, up to four pumps can be combined in one system.

The entire range from micro- to semi-preparative HPLC is covered with our assortment of different pump heads, also available as bio-compatible options.

Smartline Manager 5000

This combination instrument can be equipped with up to 3 modules: a quaternary low pressure gradient valve module, a four-channel degasser with innovative membrane technology (Teflon® AF), and an interface board for integration of analog or third-party instruments.

Smartline PDA Detector 2800

This diode array detector covers the entire wavelength range from UV through VIS up to the near IR by combining a deuterium and a tungsten/halogen lamp.

It has a very high resolution at 0.8 nm per pixel and functions with a wavelength accuracy of < 0.5 nm and a wavelength reproducibility (precision) of < 0.1 nm.

Three-dimensional data can be recorded to discover full spectral information for the peaks in your chromatograms as well as the spectra of each of the substances separated. If you only perform measurements in the UV range, this detector is also available as a one lamp device.

The Smartline PDA Detector 2800 is equipped with fiber optics, making it possible to spatially separate the measurement cell from the detector.

Smartline UV Detector 2600

Our multiple-wavelength detector based on PDA technology is particularly impressive for its outstanding noise and drift specifications. The wavelength scanning speed (0.1 sec/scan) is exceedingly fast.

If you want to acquire data for more than one wavelength per run, this is the detector of choice.

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Smartline – HPLC by Design

Smartline UV Detector 2500

This wavelength-programmable routine detector is ideal for use in micro-HPLC, analytical, and preparative HPLC with flow rates of up to 10 l/min. As with all KNAUER detectors, the Smartline UV Detector 2500 is very robust and service-friendly. The optical bench can be easily pulled out from the front of the instrument: a lamp exchange has never been this easy! An alternative model with fiber optics is available for measurements at extreme temperatures or in hazardous environments.

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Smartline RI Detector 2300 / 2400

This detector is the best choice for detecting substances not absorbing in the UV range such as carbohydrates or for the analysis of polymers in size exclusion chromatography. The analytical model of this detector (RI Detector 2300) is designed for flow rates up to 5 ml/min and the preparative model (RI Detector 2400) for flow rates up to 100 ml/min.

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Smartline Autosampler 3900

The autosampler for more sophisticated applications. Standard features include partial loopfill and precolumn derivatization capability.

The injection volume can be selected from 1–5000 µl in 1 µl increments.

The Smartline Autosampler 3900 is also available with a preparative option (3900 Prep), tray cooling (3900 Cool), a biocompatible option (3900 Bio), or combinations of these.

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Smartline Autosampler 3800

Our universal autosampler for routine analyses. It can hold up to 96 standard 1.8 ml vials in its sample carousel.

Fixed volumes from 5–5000 µl can be injected with high precision and accuracy by selecting appropriate sample loop.

This autosampler is also optionally available with partial loopfill operation (3800 Plus), sample tray cooling (3800 Cool) and as a biocompatible model (3800 Bio).

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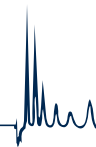
Smartline Column Oven 4000

Optimally suited for the majority of HPLC applications. This oven features heating and active Peltier cooling in the range from 10° below room temperature up to 140 °C. Its contact oven design and thermostating of the mobile phase before the column inlet ensure an even temperature distribution over the entire column for improved separation results. Active column eluate cooling minimizes peak broadening and noise/drift in subsequent detection

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Everything under Control

ChromGate® Software

ChromGate is a Client/Server structured 32-bit software and provides the highest available level of data security and GLP compliance.

It fully implements the guidelines specified by FDA 21 CFR Part 11 for the storage and protection of electronic archived data and electronic signatures.

Since the software is sold as license options, you only buy the software that you really need.

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Other KNAUER Products for Chromatography

Preparative HPLC Systems

KNAUER offers complete systems for the reliable purification of value products by HPLC. Pressure stability up to 400 bar, flow rates up to 1000 ml/min, eluent recycling, as well as a variety of detection principles are available. All components are optimally matched to each other and can also be controlled through software, enabling precise fraction collection and unique features. GLP compliance is also included.

Simulated Moving Bed Systems (SMB)

KNAUER has many years of experience with this chromatography technology for the continuous extraction of pure substances.

SMB from KNAUER is used with great success in many areas of chemistry and biochemistry. Compared to preparative HPLC, one can save up to 90% of the eluent and column packing materials are also more economical. Up to 1,000 kg of pure substance can be produced per year by using a KNAUER SMB system.

HPLC Columns

KNAUER offers separation columns from 2–62 mm inner diameter, precolumns as well as stationary phases for a large variety of applications, including high-performance phases such as our Eurospher. Every column is individually tested to ensure the highest quality and reproducibility of your results.

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Visit www.knauer.net for details on complete HPLC systems, HPLC columns, and osmometers.

KNAUER · ASI · Advanced Scientific Instruments
Wissenschaftliche Gerätebau
Dr. Ing. Herbert Knauer GmbH
Hegauer Weg 38 · D-14163 Berlin, Germany
Telephone: +49 (0)30 / 80 97 27-0
Telefax: +49 (0)30 / 8 01 50 10
E-Mail: info@knauer.net
Internet: www.knauer.net

Your local distributor: