### Science Together

## JUL KNALER

## THE INFINITE COLUMN - PEAK RECYCLING ELONGATES THE SEPARATION LENGTH OF THE COLUMN


Christian Benkhäuser, Hagen Schlicke; benkhaeuser@knauer.net KNAUER Wissenschaftliche Geräte GmbH, Hegauer Weg 38, 14163 Berlin; www.knauer.net


#### **SUMMARY**

Peak recycling basically elongates the separation length of the column by using the column bed multiple times. Most preparative chromatography users are familiar with the term but do not necessarily know in which cases this special solution is beneficial. The technology allows the eluted substances to be led back to the column after these have passed through it. Thereby, critical peak pairs that run together can be separated further with each cycle. The only requirement for peak recycling is an isocratic separation method, one additional valve, and a T-fitting.


#### INTRODUCTION

A preparative HPLC system consist of an injection system, a preparative pump, a preparative column, a detector, and a fractionation valve. Chromatography columns for preparative applications are often cost intensive. Hence, the user intends to find the best compromise between incurring costs and the obtained separation efficiency. Thereby, the individual separation challenge and the requirements for the purification set the conditions. That is why complex separation challenges can lead to high expenses and cost-ineffective solutions. In cases where special phases or small particle sizes are required for an efficient purification, not only the cost factor of the preparative column but also the system requirements are limiting. Peak recycling offers a surprisingly simple solution to increase the separation performance without the need for larger investments.

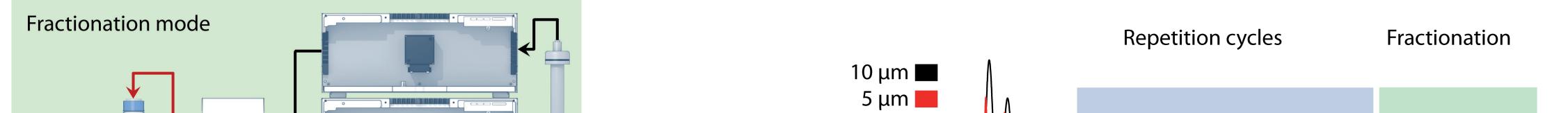

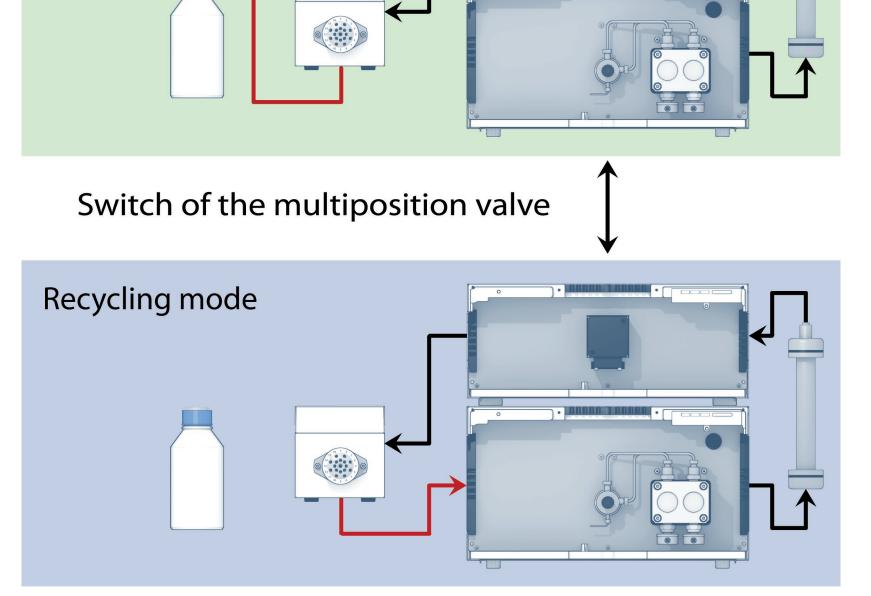

**Fig. 1** Preparative HPLC system with multiposition valve enabling peak recycling.

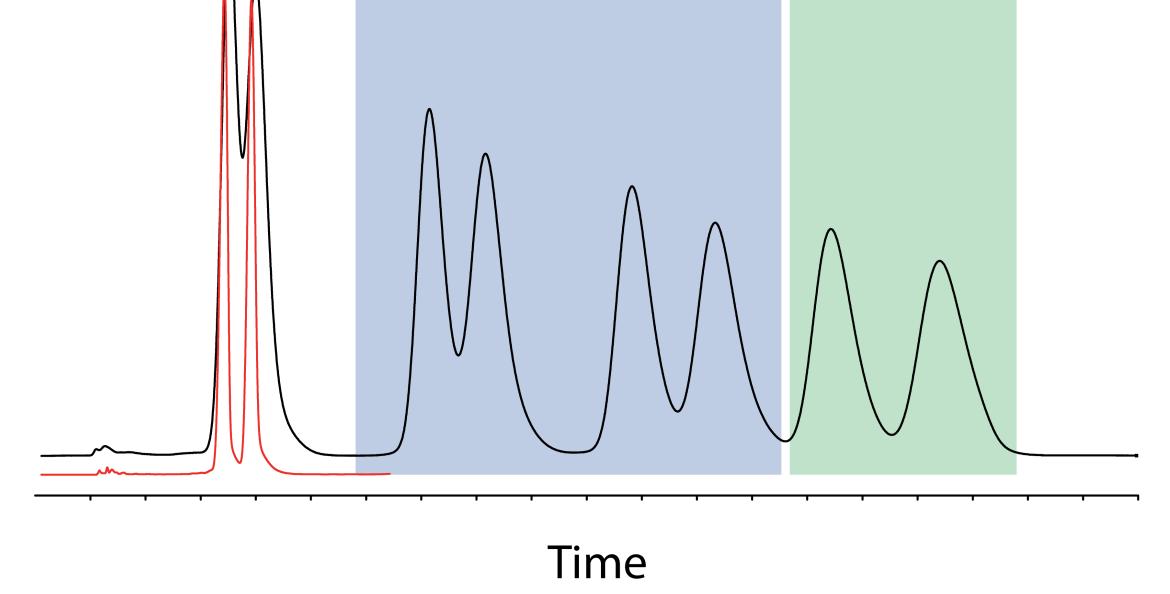

**Fig. 2** Comparison of the separations of  $(+/-)-\alpha$ -Tocopherol with more (5  $\mu$ m, red) and less expensive column material (10  $\mu$ m, black) with the help of peak recycling.

### **TECHNICAL REQUIREMENTS**

To implement the recycling mode merely a multiposition value and a T-fitting are required. The value is connected by the T-fitting to the system pump supply via a port (in most cases the second last port of the value). The value is switched to that particular port when the system is supposed to skip from normal to recycling mode. Subsequently, the pump will draw the eluent not from the stock but rather recycle the solvent through the entire system (**Fig. 1**).

The sample will be pumped in cycles until the valve is switched to fractionation or waste position.

#### **APPLICATION**

The separation of the two (+/-)- $\alpha$ -Tocopherol isomers is a practical example for the optimization of a separation procedure with the help of peak recycling. For preparative HPLC the potential for optimization is far more limited than for analytical HPLC. In reversed phase mode, the choice of column materials is practically restricted to standard phases with particle sizes of usually 10  $\mu$ m. Such preparative columns with inner diameters of 20 to 100 mm show on the one hand a lower separation performance but on the other hand cause a much lower back pressure at high flow rates, than analogous columns with small particle sizes. Users who are facing a complicated separation task like tocopherol and its isomers, where the target substance elutes narrowly to the contaminant, will be quickly limited in their options (**Fig. 2**). The costs for special phases leading to better separation and the required high performance pumps often exceed the benefit. Anyhow, a compromise of normally impaired separation performance is not even necessary. Using the the right technique, great separation results can also be achieved with such inexpensive large particle materials. For the purification of (+/-)- $\alpha$ -Tocopherol at standard conditions, usually, the user has to pay for higher yield with lower purity or vice versa. However, by the elongation of the separation length by peak recycling a baseline separation can be achieved already after three repetition cycles (**Fig. 2**, black). Fractionation can hence be carried out without compromising purity nor yield.

#### CONCLUSIONS

Peak recycling (**Fig. 2**, black) is able to enhance the separation of the (+/-)-α-Tocopherol isomers. Purity and yield are comparable to the separation with much more expensive column material (**Fig. 2**, red).

A standard HPLC system can be easily converted for the implementation of peak recycling. The modular architecture of KNAUER HPLC systems makes them ideal for the implementation and optimization of such techniques. Components like the multiposition valve for switching between the fractionation and recycling mode can be easily integrated into the system.



Additional information:

VTN0001 © KNAUER Wissenschaftliche Geräte GmbH

### Science Together

## **MKNAUER**

# THE INFINITE COLUMN - PEAK RECYCLING ELONGATES THE SEPARATION LENGTH OF THE COLUMN

#### **ADDITIONAL INFORMATION**

The terms "recycling" as well as "recycling mode" are commonly used to describe different types of recycling:

1. Time-based peak recycling:

This mode is not automated, applicable with any chromatography software, but can be more tedious to configure. Shifts in retention time within one separation can cause differences in purity of the collected fractions.

2. Automated peak recycling:

With this mode the software controls the recycling by peak height and/or slope. Particular software packages are needed for the automated mode (e.g. OpenLAB CDS EZChrom Edition) offering high level of comfort and safety.

3. Solvent recycling:

Some software packages like ClarityChrom Prep or PurityChrom offer the so-called solvent recycling as an additional option to save solvent. If the solvent in chromatogram regions without peaks can be considered clean, the solvent is recycled back into the eluent stock. This can be either based on detector signal or time. This mode can considerably lower the operating costs of a system.

#### **ADDITIONAL MATERIALS AND METHODS**

#### Tab. A1 Method parameters

Eluent	Hexane/2-Butanol (99:1, v/v)		
Gradient	isocratic		
Flow rate	20 mL/min		
Column temperature	RT	Run time	15 min
Injection volume	0.2 mL	Injection mode	Manual
Detection wavelength	290 nm	Data rate	10 Hz
		Time constant	0.1 sec

Tab. A2 System configuration & data

Instrument	Description	Article No.
Pump	AZURA P 6.1L isocratic, 50 mL pump head	APH30FA
Valve	Manual 6-port/3-channel injection valve, 1/16", SSt	A1357
Sample loop	Sample loop, 1/16", Sst	A0565
Assistant	AZURA ASM 2.1L Left: AZURA UV Detector UVD 2.1S Middle: AZURA V 2.1S with 16-port multi-position valve head Right: -	ADA00 AWA30BH
Flow cell	Semi-preparative UV Flow Cell, 1/16"	A4042
Column	Cellulose-modified silica phases for enantiomer separation (20 mm x 250 mm)	On request
Software	PurityChrom <sup>®</sup> 5	A2650

