

## Fast Separation of organic solvents

**Method** VEV0033J  
**HPLC**

**Column:** Eurokat H, 10 µm, 120 x 8 mm

**Order No.** 11GX340EKN

**Phase:** Eurokat H, 10 µm

**Conditions:** Eluent: A: 0.01 N Sulfuric acid  
Flow rate: 1.5 ml/min  
Temperature: 75 °C  
Volume: 10 µl

**Detection:** RI

**Substances:** Acetonitrile, Dimethylsulfoxide, Tetrahydrofurane, Dimethylformamide

**Keywords:** Organic solvents, Eurokat H

**Chromatogram:**

- 1 Acetonitrile
- 2 Dimethylsulfoxide
- 3 Tetrahydrofurane
- 4 Dimethylformamide

