

## Determination of Carbohydrates with Eurokat Pb

**Method** VFD0139J  
**HPLC**

**Column:** Eurokat Pb, 10 µm, 300 x 8 mm

**Order No.** 30GX350EKN

**Phase:** Eurokat Pb, 10 µm

**Eluent:** Water  
**Gradient:** Isocratic  
**Flow rate:** 0.5 ml/min  
**Temperature:** 75 °C  
**Volume:** 50 µL

**Detection:** RI

**Substances:** Glucose, Sucrose, Lactose

**Keywords:** Lactose, Eurokat Pb, Carbohydrates, Sugar

**Chromatogram:**

- 1 Sucrose
- 2 Lactose
- 3 Glucose

