

Release Notes

KNAUER Drivers 7.2.0.1038 for Chromeleon 7.2

Version 1.6, May 2017

Thank you for choosing KNAUER drivers for Chromeleon 7.2. This document describes important information about the installation and use of the KNAUER drivers and extensions for Chromeleon 7.2. Please read this document carefully before you install and start using this product.

System Configuration

Version Information

The KNAUER drivers require a Chromeleon Version 7.2. With other versions, the KNAUER drivers and extensions may not work properly.

Supported KNAUER Instruments and minimum Firmware Version

Pump P 2.1L	FW V01.07.000-P0023
Pump P 6.1L	FW V01.02.000-P0052
Detector UVD 2.1L	FW V02.04.000-P0036
Detector MWD 2.1L	FW V01.01.000
Detector DAD 2.1L	FW V01.01.000
Detector DAD 6.1L	FW V01.02.000
Detector RID 2.1L	FW V01.13.00.000
Autosampler 3950	FW V1.15
Autosampler AS 6.1L	FW V1.17
Column Thermostat CT 2.1	FW V01.06

System Requirements

Component	Minimum Requirement	Recommended Requirement
Processor	3.0 GHz Intel Core 2 Duo	3 GHz Intel Core i7 or better
RAM	3 GB	8 GB or more
HDD	60 GB	120 GB for systems with DAD Detectors
Optical Drive	DVD	
Display	1024 x 768 resolution 32-bit color	1280 x 1024 or higher 32-bit color
USB ports	1 port for license USB stick	
Ethernet port	1 port for LAN router	

OS Requirements

Microsoft Windows 8.1

Professional Edition 64-bit version

Microsoft Windows 7

SP1 Professional or Ultimate Edition, 64-bit version

Using of 32-bit operating system is not recommended

Microsoft Windows Vista

SP2 Business or Ultimate Edition, 32-bit Version

Microsoft Windows Vista is not recommended

Driver Requirements

KNAUER drivers for Chromeleon 7.2 will require a Chromeleon 7.2 Instrument Controller Option – Instrument class 3.

Release History

April 2016 KNAUER Drivers 7.2.0.1024

April 2017 KNAUER Drivers 7.2.0.1038