



KNAUER
Wissenschaftliche Geräte GmbH
Hegauer Weg 38
14163 Berlin, Germany

Phone: +49 30 809727-0

Fax: +49 30 8015010

E-Mail: info@knauer.net


Internet: www.knauer.net



Read the corresponding technical documentation for handling and safety reasons.

Mounting the pH flow cell onto the AZURA® Organizer

(KNAUER Order no. A70085)

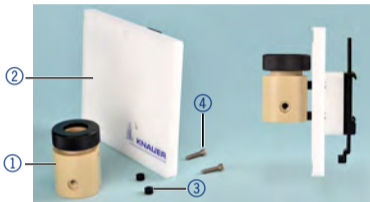

1. Remove the black holder from the pH flow cell ①.
2. Put the two screws ④ through the holes in the organizer ②.
3. On the side of the flow cell place the spacers ③ between flow cell and organizer.
4. Hand tighten the screws using the enclosed L-wrench.