




Read the corresponding technical documentation for handling and safety reasons.

Installation capillaries 1/16"

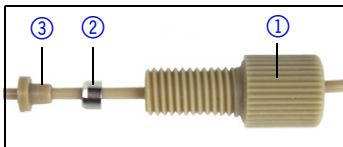

1. Ensure that the capillary ends are flat and even.
2. Push the capillary through the head of the nut ① .
3. Slide the seal ring ② over the capillary. The tapered end has to point away from the nut.
4. Attach the ferrule ③ .
5. Fasten the capillary hand-tight.



KNAUER

KNAUER

Wissenschaftliche Geräte GmbH

Hegauer Weg 38

14163 Berlin, Germany

Phone: +49 30 809727-0

Fax: +49 30 8015010

E-Mail: info@knauer.net

Internet: www.knauer.net