Azura

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Table of Contents

NOTE! For your own safety, read the manual and always observe the warnings and safety information on the device and in the manual!

Mounting the Mixer	4
Dismounting the Mixer	5
Repeat Orders	6
Technical Data	6

Mounting the Mixer

You mount the mixer with 2 Allen screws. The screws are included with the accessory kit.

Tools

- Allen screwdriver, 2 mm
- torque wrench, 2.5 mm

Hinweis

The biocompatible mixer can be used with stainless steel or PEEK fittings for the capillaries. In case you use stainless steel, the sealing ring must be made from PEEK.

CAUTION!



Damage to components caused by overtightened fittings!

Note the torque of the fittings:

- For stainless-steel fittings: 5 Nm
- For PEEK fittings: maximum 1 Nm or finger-tight and an additional quarter of a turn with appropriate pliers.

Process and Figure

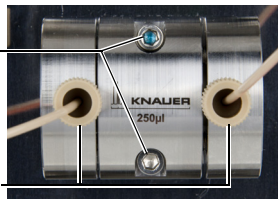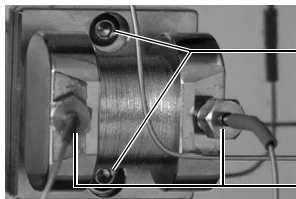1. Attach the mixer and fix it with the screws ①.
2. Using the screwdriver, fasten the screws.
3. Fasten the screw fittings ② together with the capillaries.

Afterwards take the device into operation.

Dismounting the Mixer

The mixer is dismounted in one piece, you are not allowed to disassemble the subassembly.

Prerequisites

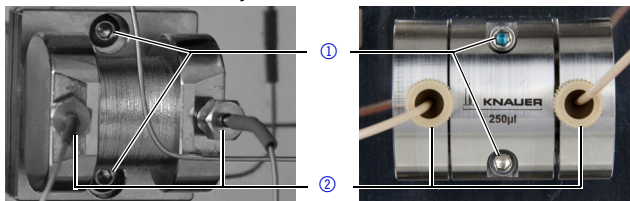
- The mixer has been flushed with ethanol.
- The hole plugs are laid out ready.

Tools

- open-end wrench, sizes 1, 4
- Allen screwdriver, 2 mm

Process


1. Using the open-end wrench, unfasten the fitting ②.
2. Close the openings with the hole plugs.
3. Using the screwdriver, unfasten the screws ①.
4. Remove the mixer and lay it to the side.

Connect the capillaries, so there is no way for solvent to leak. Afterwards take the device into operation.

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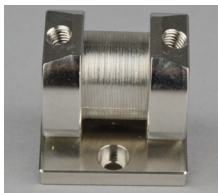KNAUER
Wissenschaftliche Geräte GmbH
All rights reserved.
The information in this document
is subject to change without prior
notice.
Translation of the original German
edition of this manual, version 1.0
Last updated: 2015-05-04
Printed in Germany on environmen-
tally friendly paper from sustainable
forests.

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KNAUER Wissenschaftliche Geräte GmbH