

Information for Installation of an Offline Version of PurityChrom® MCC Plus

1. PurityChrom MCC Plus operates under Microsoft Windows XP, Windows Vista, Windows 7, Windows 8, and Windows 10.
2. The user account settings have to be set to “Never notify” before installation. After the installation, the settings can be reset. (Start → Control Panel → User Accounts → User Accounts → Change User Account Control Settings → Never notify)
3. Restart your computer.
4. Insert the CD which includes PurityChrom MCC Plus and the configuration files of your system.
5. Start the installation as an administrator. The installer will ask you to choose the language of the software, your integrated UV-detector, and the destination location for the files that will be installed. Do not change the given destination. Choosing “Demo-Version” in the installer is not necessary because when PurityChrom MCC Plus will be opened without license file and dongle, the software switches automatically into demo mode.
6. If the installation is finished, copy and replace the PurityChrom.ini-file from the computer of your active FPLC system (located under *C:\Windows*) or from the PurityChrom MCC Plus CD (folder: *INI*) to *C:\Windows*.
7. Copy and replace the PurityChrom.cfg-file from the computer of your active FPLC system (located under *C:\PurityChrom-MCC-Plus*) or from the PurityChrom MCC Plus CD (folder: *CFG*) to *C:\PurityChrom-MCC-Plus*.
8. Copy the background (*.bmp) from the computer of your active FPLC system (located under *C:\PurityChrom-MCC-Plus\Visualisation\Visualisation Backgrounds*) or from the PurityChrom MCC Plus CD (folder: *Visualisation\Background*) to *C:\PurityChrom-MCC-Plus\Visualisation\Visualisation Backgrounds*
9. Copy the visualisation (*.vis) from the computer of your active FPLC system (located under *C:\PurityChrom-MCC-Plus\Visualisation\Visualisation Files*) or from the PurityChrom MCC Plus located on the CD (folder: *Visualisation\Files*) to *C:\PurityChrom-MCC-Plus\Visualisation\Visualisation Files*
10. **Installation under Windows versions older than 8 and 10:** Install PurityChrom MCC Plus. Go to *C:\PurityChrom-MCC-Plus* and right-click the PurityChrom-MCC-Plus.exe to open the properties. Go to compatibility and check “Run this program in compatibility mode for” and choose “Windows Vista (Service Pack 2). After this, check “Run this program as an administrator“ click ”Apply“ and ”Ok“, then close the window.
11. **Installation under Windows 8 or 10:** Install PurityChrom MCC Plus. Go to *C:\PurityChrom-MCC-Plus* and right-click the PurityChrom-MCC-Plus.exe to run the program as administrator. Open the visualization and the communication log once.

12. To open result or method files from your existing system, simply transfer them to *C:\PurityChrom-MCC-Plus\Data*.
13. If you use the user administration of PurityChrom MCC Plus and want to have the same settings in your offline version, then copy and replace the file *UserAdminsitration.cfg* from the computer of your active FPLC system (located under *C:\PurityChrom-MCC-Plus*) to *C:\PurityChrom-MCC-Plus*.